



1 WHAT'S IN THE BOX

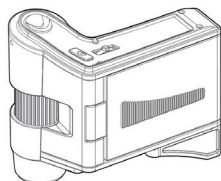

- FlipView® LCD digital handheld microscope
- Full metal stand with 1/4-20 thread
- AC adapter
- Lithium ion battery (1050 mAh)
- USB cable
- Video out cable for connecting to external monitor
- Quick start guide
- Screen/lens cleaning cloth
- Calibration ruler
- Carry bag


2 INSTALL THE BATTERY

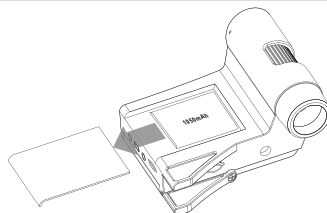

1. Open battery door by sliding the door outwards.
2. Insert lithium ion battery.
3. Make sure the batteries' polarities are correct.

NOTE: Always fully charge your 1050 mAh lithium battery before using.


3 INSERT YOUR MICRO SD CARD

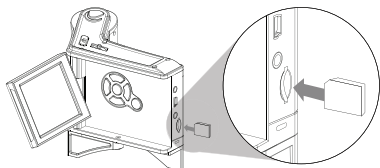

1. Insert Micro SD card into Micro SD card slot.
2. Press gently until the card clicks into place.


NOTE: Micro SD card not included. You can view specimens without a Micro SD card, but cannot record video or save images.

4 USING YOUR MICROSCOPE

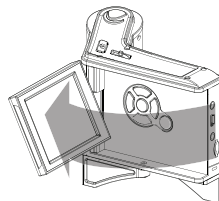


1. Flip open the LCD screen.
You are now ready to enjoy your FlipView LCD Handheld microscope!
NOTE: If the screen turns off while open due to energy save setting, press the ON/OFF button to power ON.
2. When you are finished viewing, flip the LCD screen closed to turn off.

5 CONNECT TO YOUR COMPUTER

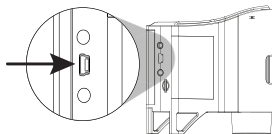

1. Plug the USB cable into an available USB port on the computer (Windows or Mac).
2. Plug the other end of the same cable into the FlipView mini USB port.
3. The LCD screen on the FlipView will display 2 options:
A. Upload Images (transferring images to PC)
B. Preview on PC (using included software)
4. Make your choice and refer to the downloadable manual for detailed instructions.

6 CONNECT TO YOUR TV

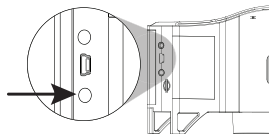

1. Insert the 1/8" plug end of the video out cable into the AV out port on FlipView.
2. Insert the RCA plug on the video out cable into the TV/monitor's video in ports.

NOTE: The video in ports on most displays and televisions are color coded yellow.

NOTE: Download and install software for the PC from the product page at celestron.com.